Fiche technique

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Vertical shower (ceiling mounted) – ref 1100

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This ceiling mounted shower minimizes required space with a water inlet by ceiling.
Pull-rod activated.

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Material

Galvanized steel pipe.
Covered by a PL11 polyamid resin (anti-corrosion),
Stainless Steel valve and brass connections,
ABS Shower head (Option Stainless Steel)
Pull-Rod steel
Or All stainless steel model 304 to 316.

Water inlet/output

Water inlet 1" BSP
Pressure : 2 bars recommended - Min 1.5 Max 8

Shower

Manual activation by pull rod.
(anti-kickback system)
Flow: 110L/min for a pressure between 2 and 8 bars.
Valve 1" with pressure regulation

Plate

Shower plate
Photo luminescent with green background
AFNOR X08-003 Standard

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Normes / Standards
EN 15154-1, DIN 12899-3:2007,
ANSI/SEA Z358: 2009

Options

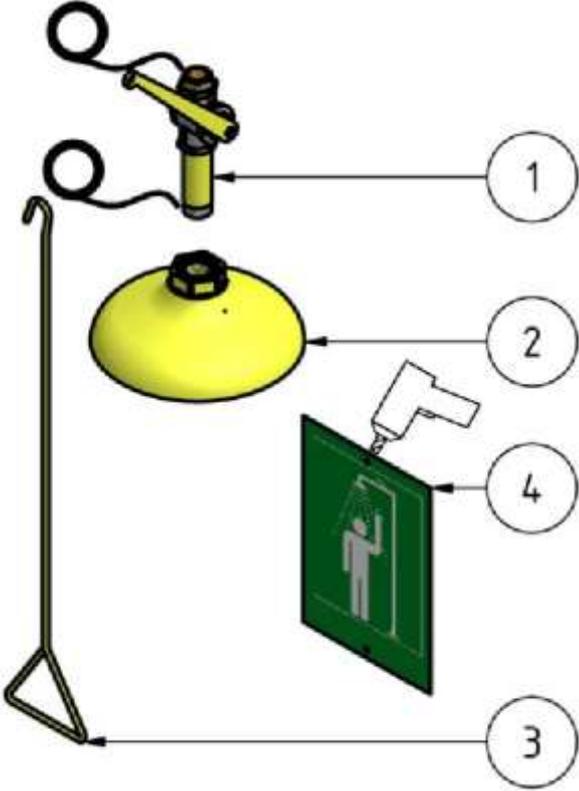
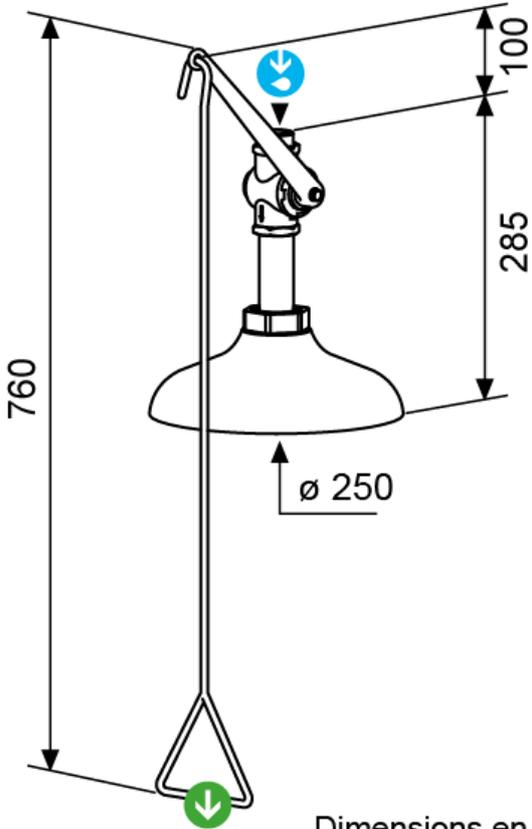
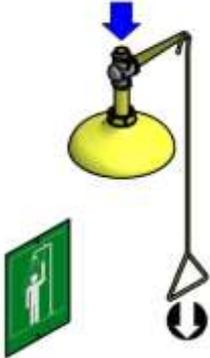
Pomme inox *Stainless steel shower head*

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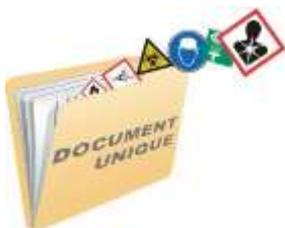
*Install these security equipment in order to make them **visible and accessible from at least 3 ways**.
The emergency equipment must be easily accessible by the user (**less than 10 sec**).
A correct operation test must be planned **every week** by an internal procedure.*

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Installation and use for showers and eyewashes. Devices Connections

1. The emergency decontamination booth must be close to danger exposure. **(Max 10 sec or 30M obstacle)**.
*In case of high probability of simultaneous accidents, **anticipate more than only one equipment**.*
2. Safety equipment must be **easily accessed and clearly identified**
3. Safety Equipment must be connected to the drinking water system with stop valve for maintenance.
4. Test must be executed **every week by an internal procedure**.
5. For an outdoor facility, use frost-protected or freeze-proof equipment.
6. All the personnel must know how to use correctly security equipment.
7. **Eyewash time: from 10 to 15 minutes.**
8. **Shower time: from 5 to 10 minutes.**

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